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REV. 1A12M11

INDEX

Keyboard and keys function	4
Power Off	4
Zero and Tare functions	4
Unit weight selection	4
One point calibration	5
Two point calibration	5
Auto-power off function	6
Division selection	6
High and low limits setting	6
Error messages	7
Battery instructions	8
Guarantee	8
EC Declaration of Conformity	20

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KEYBOARD

KEY FUNCTION

Switch on / off the scale.
Zero and Tare key.

To select unit weight
To confirm in calibrate procedure.

Switch function

Position choose function

OPERATION

Power Off

Press the  key for 2 seconds, after display “SLEEP”, then release the key to power off the scale.

Zero and Tare

Press the  key one time for zero function, zero range is 4% of full capacity.
If the weight is over this range, it will make the Tare operation, and the tare indicator will light on.

Unit weight selection

Press the  key to select the desired unit weight (lb,oz,g)

CALIBRATION

One point calibration.

1. Under normal weighing mode, press  for 2 seconds until display shows "SLEEP".
2. Press the  key to catch zero, the display will show "-----" and keep the tray stable.
3. Display: xxx, clue to load xxx g poise, now you can press  to choose the load point, you can load on the poise which you have choose, press the  key to confirm and keep it stable.
4. If the weight is not correct after calibration, you can try two point calibration.
Note: The two calibration methods are based on unit gram.

Two point calibration.

1. Enter to calibration mode, press the  key for 2 seconds until display shows "SLEEP"
2. Press the  key again to enter into calibration mode.
3. The display will show "F-XX" to choose capacity, press the  key to choose and press the  key to confirm.
4. Display will show: d=x to choose division, press the  enter division and key  confirm.
5. Display will show: CAL- to catch zero, empty the tray and keep it stable, then press  key to confirm, display will view "-----" catching the zero.
6. Display will show: xxx, clue to load xxx g poise, this is one point calibration, after load the poise, press  key to confirm, display will view "-----".

7. Display will show: yyy, clue to load xxx g poise, this is the two point calibration, after load the po  press key to confirm, display will view "-----"
8. If the calibration procedure is correct, the scale will return automatically to normal weighing mode.
9. If the loaded poise is low, it will display Err-3 and return to the normal weighing mode without save the calibration data.
10. During the calibration process, you can press the  to last step.

FUNCTION SETTING

Under normal weighing mode, press the  for 2 seconds, press the  to choose menu and press  key to enter in selected menu.

AUTO-POWER OFF FUNCTION

1. Display shows "S-OFF"
2. Press the  key to enter into.
3. Display shows "A-X"
4. Press the  to select the desired mode:
 - 0 = Auto-power off disable.
 - 1 = Balance will auto-power off after 1 minute of non use.
 - 3 = Balance will auto-power off after 3 minutes of non use.
 - 5 = Balance will auto-power off after 5 minutes of non use.
 - 10 = Balance will auto-power off after 10 minutes of non use.
 - 15 = Balance will auto-power off after 15 minutes of non use.
 - 30 = Balance will auto-power off after 30 minutes of non use.
 - And so on.
5. Press the  key to confirm and come back to normal weighing mode.

DIVISION SELECTION

1. Display shows “S-d”
2. Press the  key to enter into.
3. Display shows “d=xx”
4. Press the  key to choose division
5. Press the  key to confirm and come back to normal weighing mode.

HIGH AND LOW LIMIT SETTING

a) Setting alarm type

1. Display shows “S-Alr”
2. Press the  key to enter into.
3. Select the desired mode:
Alr-0= means alarm forbidden
Alr-1= means alarm in the range
Alr-2= means alarm out the range
4. Press the  key to confirm and enter to High & Low setting automatically.

b) High and Low setting

1. Display shows “ S-A-H” (High limit)
2. Press the  key to enter into.
3. Display shows “H00000”

4. Press the keys   to reduce the value.
5. Press the  key to confirm and save.
6. Display shows “ S-A-L ” (Low limit)
7. Press the  key to enter into.
8. Display shows “L00000”
9. Press the keys   to reduce the value.
10. Press the  key to confirm and save.

c) Buzzer choose.

1. Display shows “S-buz”
2. Press the  key to enter into.

d) Alarm mode setting.

1. Display shows “ S-A-S”
2. Press the  key to enter into.
3. Select the desired mode:

A-STAB= alarm when the weight is stable.
A-NSTA= alarm without stable condition.

To exit from setting mode, press the  key.

ERROR MESSAGES

- Err2, means initial zero point is not stable.
- Err3, means the loading is not enough during two point calibration.

- Err4, means the adjustment over modify range during one point calibration. Modify the range is between 0.6 times and 1.4 times of the factory calibration weight.

BATTERY INSTRUCTIONS

The scale uses AAA battery as power.

If the voltage of the batteries is low, the battery signal will appear at the left corner of display, clue to change the batteries.

The working current is 5mA, suggest to use the LR03 to work for long time.

GUARANTEE

This scale is guaranteed for one year from the delivery date. Te guarantee covers any fabrication defect of the material.

During this period GRAM PRECISION, covers the manpower and the spare parts necessary for the reparation of the scale.

This guarantee does not cover the failures caused by an inappropriate use or overcharge.

The guarantee does not cover the freight cost (transport) necessary to repair the scale.

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A large, stylized, grey 'E' logo is centered on the page. The 'E' is composed of thick, rounded strokes. The middle horizontal bar of the 'E' is a solid grey rectangle, while the vertical strokes are open at the top and bottom.

**EC DECLARATION OF CONFORMITY
DECLARACION CE de CONFORMIDAD
DECLARATION CE de CONFORMITE**

MANUFACTURER / FABRICANTE / FABRICANT:

GRAM PRECISION, S.L

ADDRESS / DIRECCIÓN / ADRESSE:

Travesia Industrial 11 / 08907 Hospitalet de Llobregat (Barcelona) – SPAIN -

PRODUCT: Electronical scale model D1

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The balances above mentioned meet with following European Standards:

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Electromagnetic compatibility 2004/108/CE

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